

Miele Standard-Wash-Programs

These programs replace the Miele-Standard-Wash-Programs described in the Programming Manual.

Program names	Wash temperature in the program blocks [°C]				Drying parameters		
	Cleaning	Sanitizing	Neutralize	Final rinse	Temperature 1 / 2 [°C]	Time 1 / 2 [min]	Cooling down [min]
INORGANICA	75		10	70	110/-	25/-	0
ORGANICA	85		10	80	110/-	25/-	0
STANDARD	70		10	70	110/ -	25/ -	0
UNIVERSAL	75		10	75	110/ -	25/ -	0
INTENSIVE	80		10	75	110/ -	25/ -	0
PLASTIC	55		10	55	70/60	10/30	0
VARIO TD	45	93 ¹⁾			90/ -	30/ -	0
SPECIAL		93	10	70	100/ -	30/ -	0
SANITIZING	85	85 ¹⁾	10		90/ -	30/ -	0
MINI	60		10	60	100/ -	30/ -	0
DI-RINSE					0/ -	0/ -	0
DRYING					90/ -	15/ -	0

1) Disinfection takes place with the final rinse.